

## Software Update notes: Blue Pippin version 6.31 / 6.40

Current software version: Linux 6.31 / Windows 6.40

### Blue Pippin

Release notes:

Date: 01/16/19

Version: 6.31 / 6.40 CD 33

#### What's new from 6.31 / 6.40 CD 31 to 6.31 / 6.40 CD 33

- Added New High Pass Plus >20kb Cassette Definition
- 

Release notes:

Date: 08/07/18

Version: 6.31 / 6.40

#### What's new from v.6.30 to 6.31 / 6.40 CD 31

- Update works for both Windows and Linux based instruments
  - Added New High Pass Plus >15kb Cassette Definition
- 

Release notes:

Date: 06/01/17

Version: 6.31

#### What's new from v.6.30 to 6.31?

- Progress Bar Update
  - Adjusted continuity test parameters
- 

Release notes:

Date: 03/09/17

Version: 6.30

#### What's new from v.6.23 to 6.30?

This update is for BluePippin users only. Under heavy pulsed-field use with BluePippin, some older electrode arrays can form a salt bridge which may interfere with performance. This update runs a diagnostic routine to check for this. The routine is run automatically at the end of every run. The routine is not visualized by the user, but a warning screen will inform users if an error is detected.

---